

# Precise Measurement of Viral Copy Number at Single-Cell Resolution for Cell and Gene Therapy Development

**Yue Wang**, Saurabh Parikh, Khushali Patel, Yilong Yang, Daniel Mendoza, Gukhan Kim, Mahir Mohiuddin\*, Hua-Jun He\*, John Elliott\*, Matt Cato, Shu Wang, Benjamin Schroeder, Chieh-Yuan Li  
Mission Bio. 400 E Jamie Ct, Suite 101, South San Francisco, CA 94080; \*National Institute of Standards and Technology. Gaithersburg, MD

## Introduction

Cell and gene therapies are transformative solutions for a host of inherited and acquired diseases for which existing interventions are ineffective. Many such therapies rely on the introduction of transgenes into host cells using viral or non-viral vectors. The accurate measurement of gene transfer is critical to the development of therapeutic agents and is a key attribute for assessing their safety and efficacy. Yet, conventional methods for measuring gene transfer lack the resolution and representation to truly reflect sample composition and either report a population average (bulk) or involve laborious and time-consuming clonal outgrowth which can take weeks. Mission Bio has developed an end-to-end solution from panel design to data analysis for single-cell targeted DNA sequencing. Here, using the Tapestri platform we demonstrate that single-cell DNA sequencing identifies transduced versus non-transduced cells with exceptional accuracy and precision, as well as measures the single-cell level vector copy number (VCN) for populations of thousands of cells with single nucleotide resolution while reducing sample processing time from weeks to days.

## Objective

To demonstrate the quantitative characterization capability of single-cell level vector copy number and transduction efficiency using the Tapestri single-cell DNA sequencing platform in terms of assay specificity, sensitivity, accuracy and precision.

## Methods

The Tapestri precision genomic platform is enabled by its novel two-step microfluidic workflow where thousands of cells are each encapsulated and lysed in the first droplet followed by cell-lysate barcoding and target DNA amplification using highly multiplexed PCR, in the second droplet. The droplets are then broken, and the amplified products are pooled for DNA library generation using Mission Bio consumables and sequenced on the Illumina NGS system. The final data is analyzed and visualized using the Tapestri Pipeline and Tapestri Insights software.

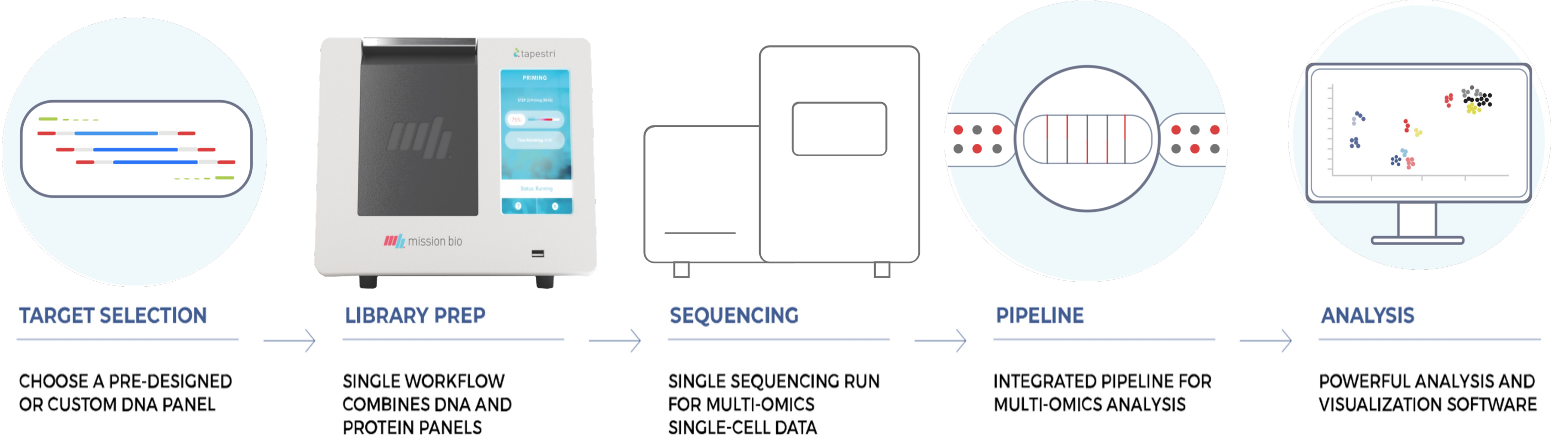


Figure 1: Tapestri workflow

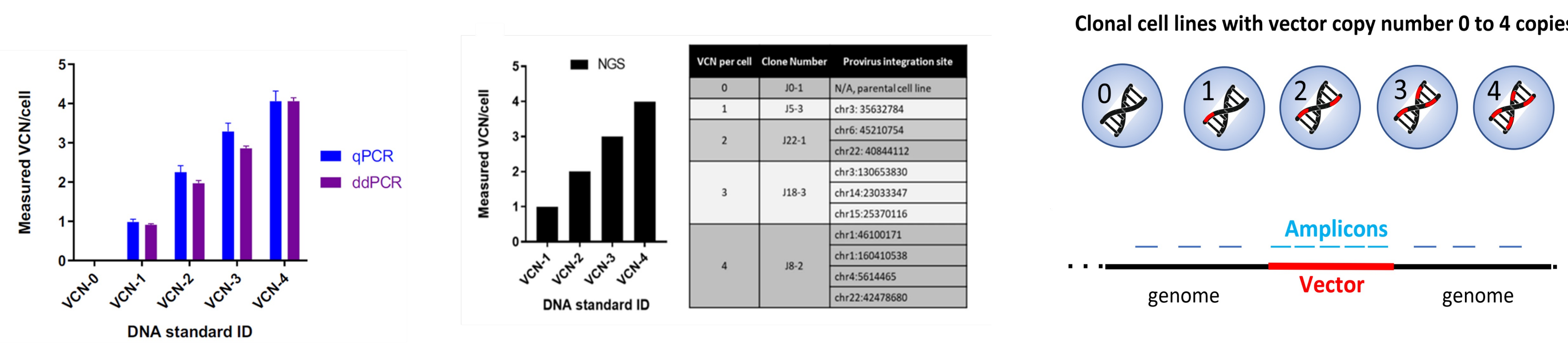


Figure 2: Lentiviral transduced VCN cell lines and vector assay design<sup>1</sup>

Characterized, experimental, lentiviral transduced Jurkat clonal cell lines with vector copy number of 0, 1, 2, 3 and 4 copies were used in this study <sup>1</sup>. A custom panel of amplicons was designed to enable the vector copy number analysis. Pure clonal cell lines of VCN 0-4 and three (3) mixtures with non-transduced Raji cell spike-in was processed for single-cell sequencing using Mission Bio’s Tapestri platform in triplicates. Tapestri data from a control cell line with known VCN, as well as unknown samples, were combined and analyzed with a copy number caller to produce single-cell level VCN calls, as well as a population average VCN and a percent transduction for each sample. A median of 5,116 cells were analyzed for each Tapestri sample.

## Results

### Tapestri vector copy number assessment tool is precise and accurate

Average copy number for various combinations of controls and test samples	Test Sample											
	VCN rep1	VCN rep2	VCN rep3	VCN rep1	VCN rep2	VCN rep3	VCN rep1	VCN rep2	VCN rep3	VCN rep1	VCN rep2	VCN rep3
Control Sample	VCN1 rep1	1.00	1.00	1.00	2.10	2.00	2.00	3.00	3.00	3.26	4.34	4.46
	VCN1 rep2	1.00	1.00	1.00	2.00	2.00	2.00	3.00	3.00	4.12	4.23	4.01
	VCN1 rep3	1.00	1.00	1.00	2.01	2.00	2.00	3.00	3.00	3.05	4.09	4.32
	VCN2 rep1	1.01	1.03	1.03	2.00	2.00	2.00	3.00	3.00	4.08	4.27	4.19
	VCN2 rep2	1.02	1.03	1.03	2.01	2.00	2.00	3.00	3.00	3.05	4.18	4.35
	VCN2 rep3	1.01	1.01	1.02	2.00	2.00	2.00	3.00	3.00	4.09	4.38	4.34
	VCN3 rep1	1.09	1.08	1.10	2.10	2.08	2.07	3.00	3.11	3.13	4.37	4.54
	VCN3 rep2	1.03	1.03	1.04	2.00	2.00	2.00	3.00	3.00	4.18	4.43	4.27
	VCN3 rep3	1.06	1.06	1.06	2.02	2.01	2.00	3.00	3.00	4.22	4.42	4.29
	VCN4 rep1	1.00	1.01	1.01	2.00	2.00	2.00	2.99	3.00	4.01	4.34	4.11
	VCN4 rep2	1.00	1.00	1.00	1.94	1.98	1.98	2.77	2.98	2.83	3.95	4.04
	VCN4 rep3	1.00	1.00	1.01	2.00	2.00	2.92	3.00	2.96	4.00	4.31	4.03

Table 1: Combinations of controls and test samples’ average VCN

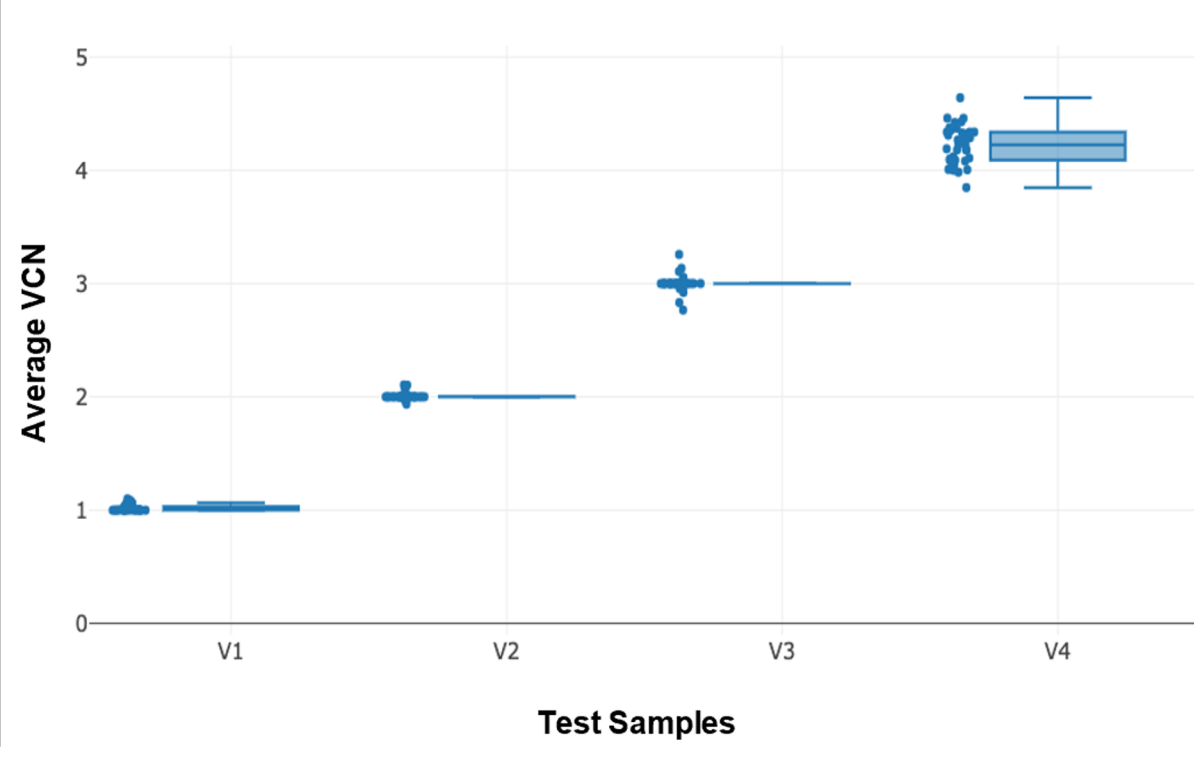


Figure 3: Control vs. average VCN

Clonal populations of engineered VCN cells (VCN1-4) were single-cell sequenced and processed through the Tapestri VCN caller individually (n=3). We demonstrate the single-cell VCN caller can use any cell population with a known number of VCNs to characterize a test sample’s VCN with high accuracy and precision (Table 1). Figure 3 is a graphical representation of Table 1, where each dot represents the average VCN of the test sample (x-axis) using various control samples. Using VCN caller model established from VCN 1 to VCN 4 each as reference points for determining average VCN produces coefficient of variation of 2.76%, 1.62%, 2.48% and 4.04%; and percent error of 2.17%, 0.43%, 0.05% and 5.56% respectively across all combinations across runs.

### Tapestri VCN measurements correlate with orthogonal ddPCR technology

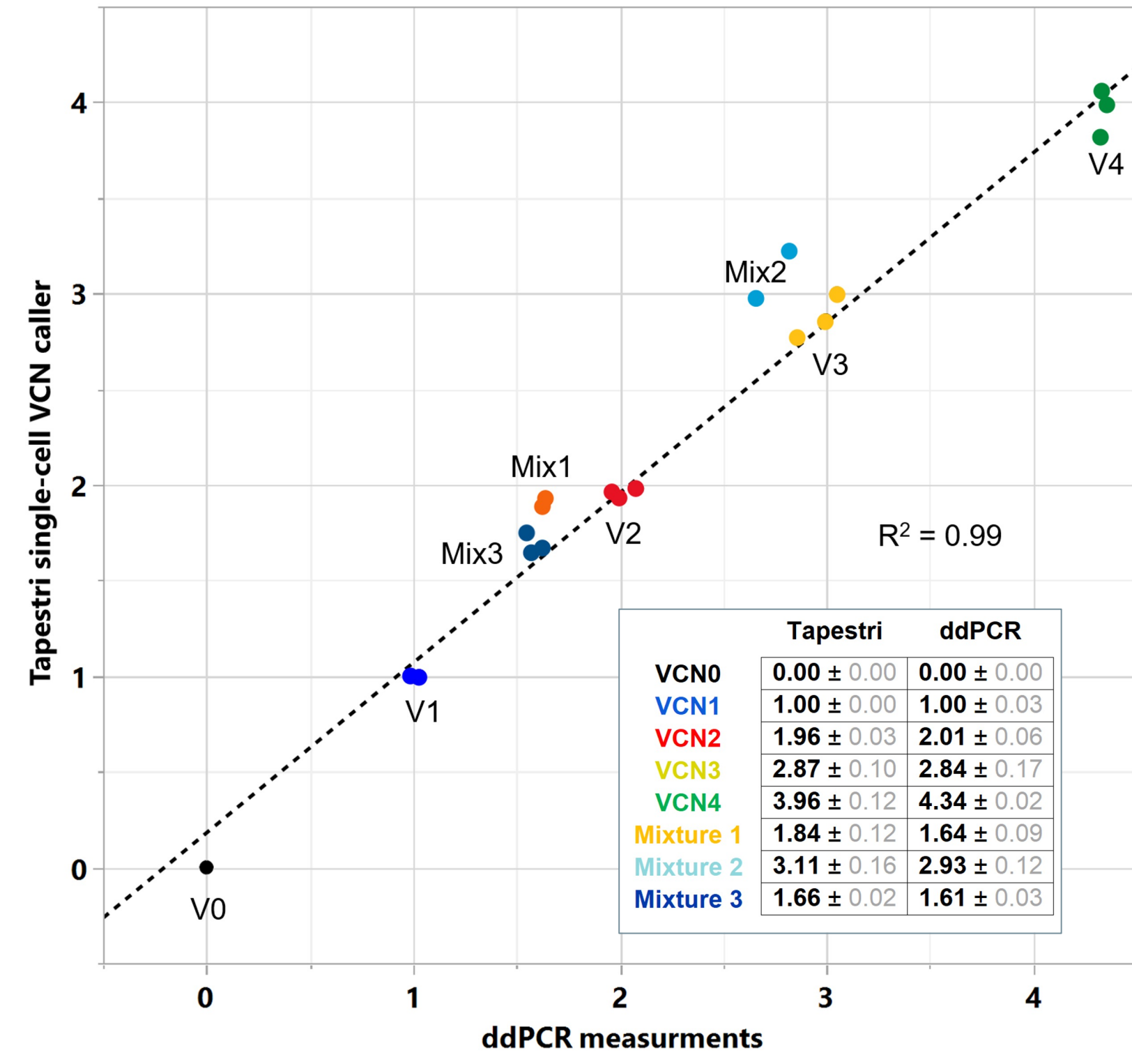


Figure 4: Tapestri vs. ddPCR VCN measurements

X-Y scatter plot of the average VCN for each sample called using Tapestri VCN caller vs. ddPCR measurement. Pure cell lines (VCN 0-4) or mixtures of cell lines (Mixes 1-3) were analyzed in triplicate by both Tapestri and ddPCR. The sample’s average VCN was then calculate based on single-cell level VCN (y-axis) and plotted against orthogonal ddPCR measurements (Bio Rad QX200 <sup>1</sup>, x-axis, n=3). The average copy number of the population estimated using Tapestri correlates well with ddPCR measurements (R<sup>2</sup>= 0.99). The average and standard deviation of replicate measurements (n=3) are shown in the bottom right table.

### Transduction efficiency can be detected with high sensitivity and specificity

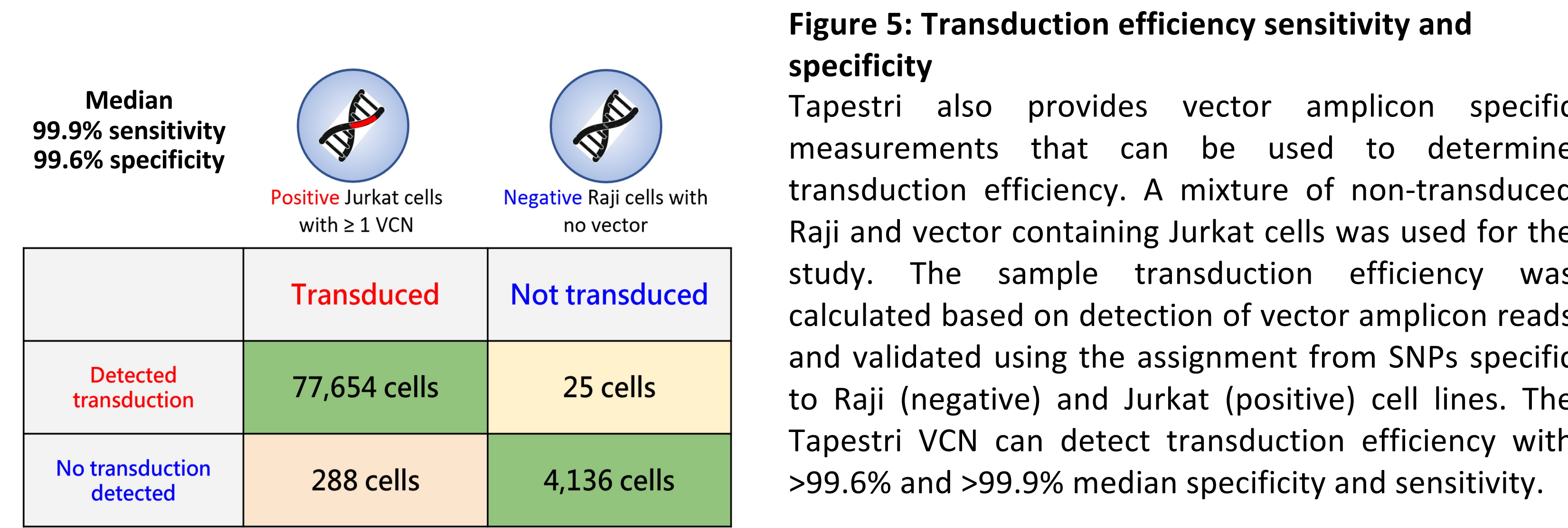


Figure 5: Transduction efficiency sensitivity and specificity

Tapestri also provides vector amplicon specific measurements that can be used to determine transduction efficiency. A mixture of non-transduced Raji and vector containing Jurkat cells was used for the study. The sample transduction efficiency was calculated based on detection of vector amplicon reads and validated using the assignment from SNPs specific to Raji (negative) and Jurkat (positive) cell lines. The Tapestri VCN can detect transduction efficiency with >99.6% and >99.9% median specificity and sensitivity.

### Single-cell level resolution of vector copy number distribution characterization

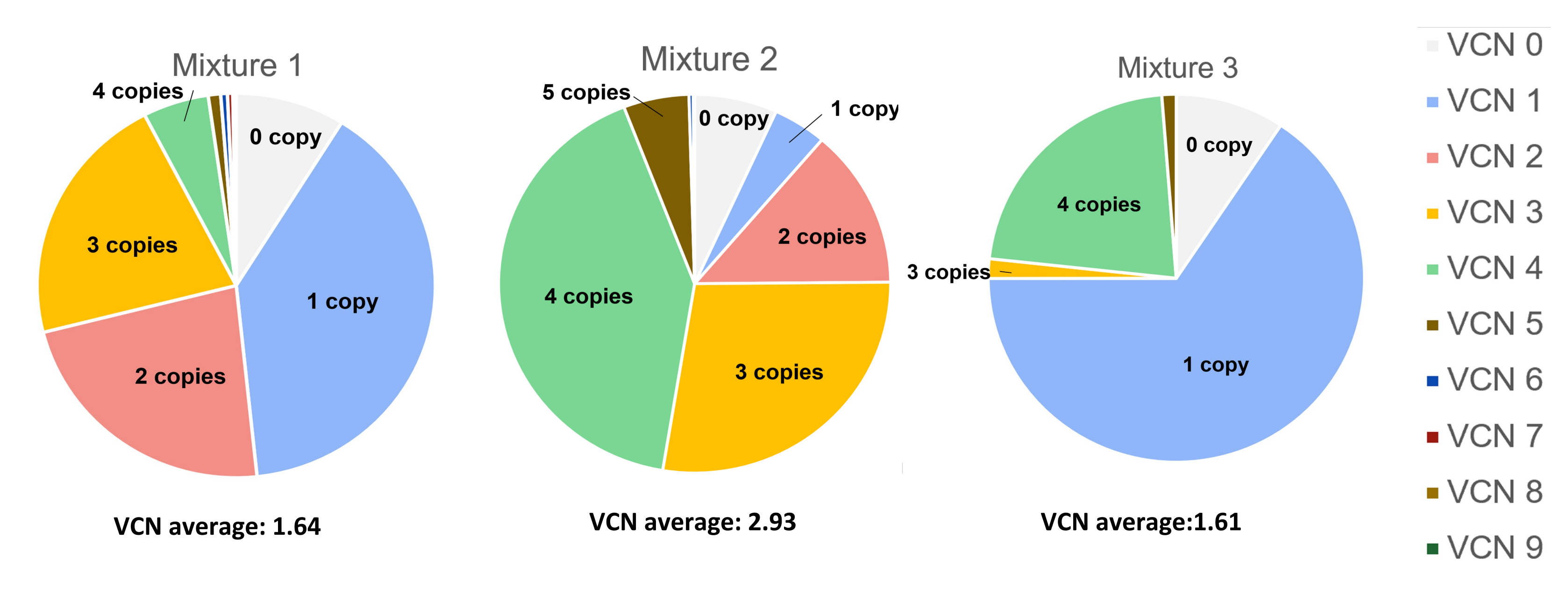


Figure 6: Single-cell VCN distribution per sample

Besides the measurement of average VCN per sample, Tapestri VCN caller calculates VCN distribution of a particular sample at single-cell level. The pie charts shows the VCN distribution calculated for each of the three mixtures, Mix1, Mix2, and Mix3 each with ~10% non-vector Raji cell spiked-in. Note that “Mixture 1” (VCN 1-4’s respective proportions = 39.4%, 22.8%, 21.3%, 5.4%) and “Mixture 3” (VCN 1-4’s respective proportions = 65.6%, 0%, 1.7%, 22.1%) VCN distributions are drastically different despite having similar average populational VCN value of 1.64 and 1.61.

### Single-cell genotype, SNPs, and transgene vector characterization

Tapestri amplicon based single-cell DNA sequencing provides the capability to analyze the genomic landscapes of regions of interest at single base resolution. The heat-map to the right shows genotyping of cells (Raji vs. Jurkat) based on unique SNPs using human amplicons. The VCN caller determines the VCN 1 vs. VCN 4 based on the read out from vector amplicons while also providing sequencing data on each amplicon. Here, Tapestri identifies a spontaneous mutation occurring in one out of four copies of the transgene vector at position 408:G/A in the VCN 4 clonal cell line, whereas VCN 1 cells contain only the expected vector sequence. Tapestri not only determines the vector copy number, but also provides sequence integrity information of transduced vector, which directly correlates with the functionality of the transgene.

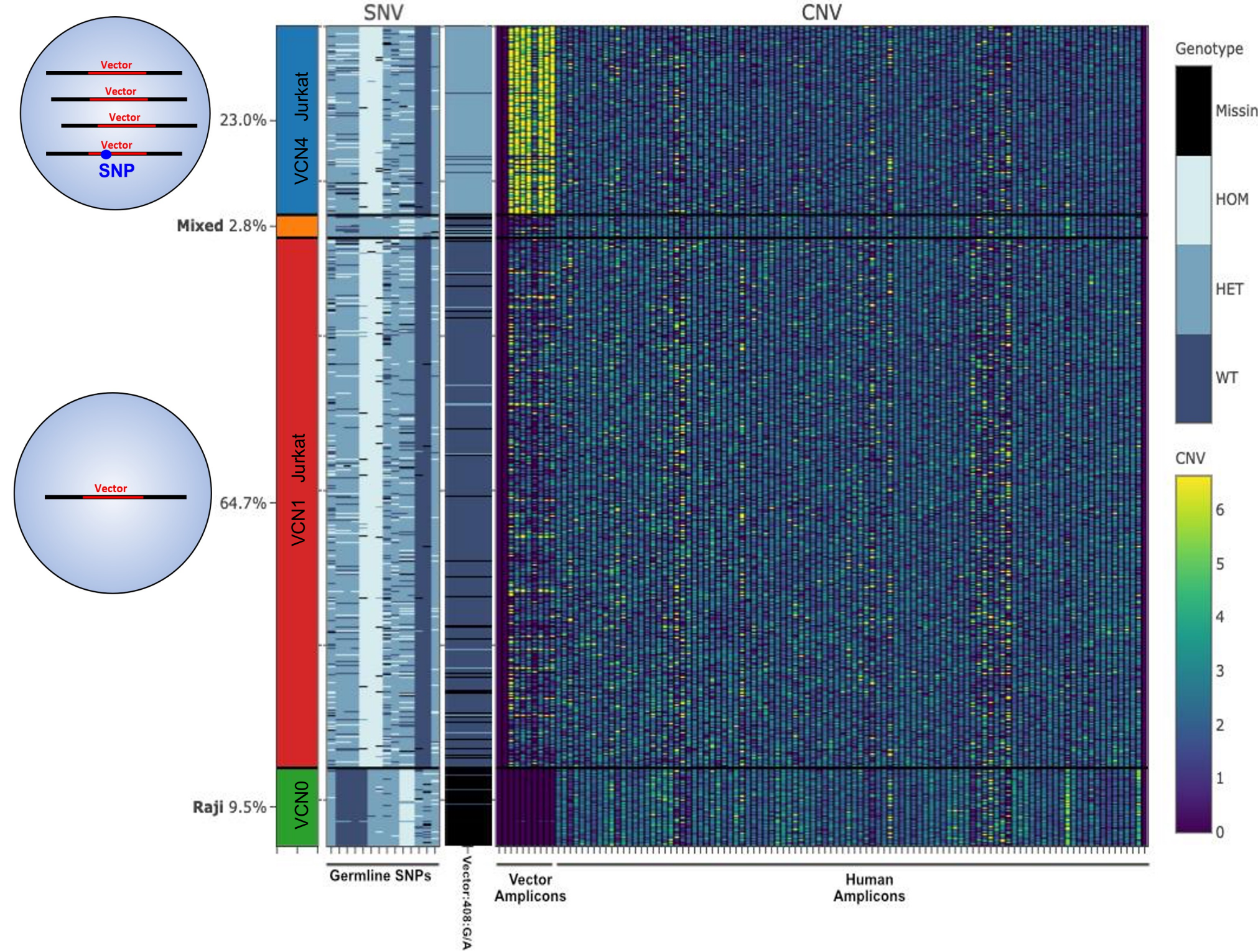


Figure 9: Heat-map of genotype, VCN, and vector sequence

## Conclusion

Using the Mission Bio Tapestri single-cell DNA sequencing platform, this study shows a consistent and reliable assay for in-depth quantification of cell and gene therapy transduction efficiency, single-cell vector copy number distribution, and populational average. The single-cell level VCN distribution, as well as single-cell vector sequence validation provides unprecedented resolution and insight to assess the potential functional efficacy and safety for CGT products. Together, these characteristics and data demonstrate the potential to accelerate and streamline both the development and release testing of cell and gene therapy products.

Reference 1. <https://doi.org/10.1038/s41598-020-79698-w>